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DirectAmp™ lysis solution

Fast Nucleic Acid Extraction from Swab, UTM, VTM and Others

DirectAmp™ lysis solution was developed for the fast sample preparation for bacterial and viral nucleic acid amplification as used in Real-Time PCR (RT-PCR), Quantitative Reverse Transcription PCR (qRT-PCR) and Colorimetric Reverse Transcription Loop-Mediated Isothermal Amplification (Colorimetric RT-LAMP).

DirectAmp™ lysis solution has been used to prepare a bacterial and viral nucleic acid sample from a wide variety of different sources, including swab, UTM, VTM and cell culture media. Using only simple heating step for lyse a sample and no nucleic acid cleanup step required towards lysis and detection.

DirectAmp™ lysis solution can be applied on a point-of-care testing to be used for massive screening that can help to isolate infectious people from the community.

Key Features and Benefits

- Simple and fast protocol reduces turnaround time (lysis as short as 10 mins)
- Minimal hands-on time and consumable required
- Affordable method with low price per sample
- Enzyme-free nucleic acid extraction
- Ideal for on-site nucleic acid testing

Applicable Specimen Types

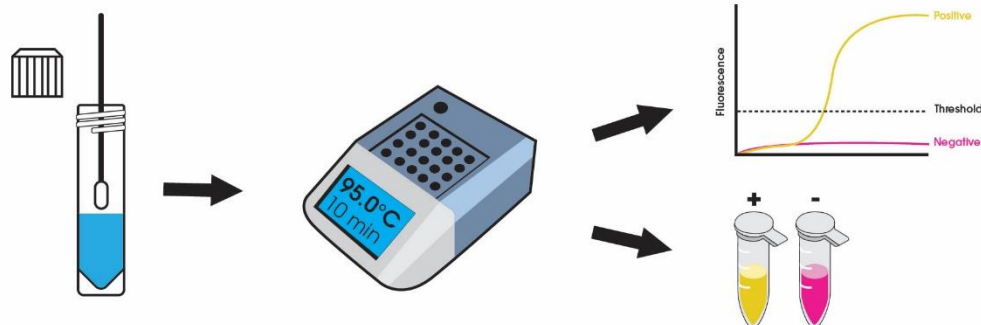
- Nasopharyngeal (NP) swab
- Oropharyngeal (throat) swab
- Nasal swab
- Oral cavity swab

Compatible instrumentation

Heating machine

- Heating block

Protocol

qRT-PCR Compatibility for COVID-19 Detection

DirectAmp™ lysis solution is compatible with nucleic acid amplification methods and does not impact on assay sensitivity or reduce fluorescence signal when used at working concentration in qRT-PCR.

Compatibility has been verified with the following kits:

- (A) Novel Coronavirus (2019-nCoV) Nucleic Acid Diagnostic Kit (Sansure Biotech, catalog S3102ESC2)
- (B) ANDIS FAST SARS-CoV-2 RT-qPCR Detection Kit (3DMed, catalog 3103010069)
- (C) Zybio Molaccu COVID-19 Detection Kit (Zybio, catalog ZB-4080102)

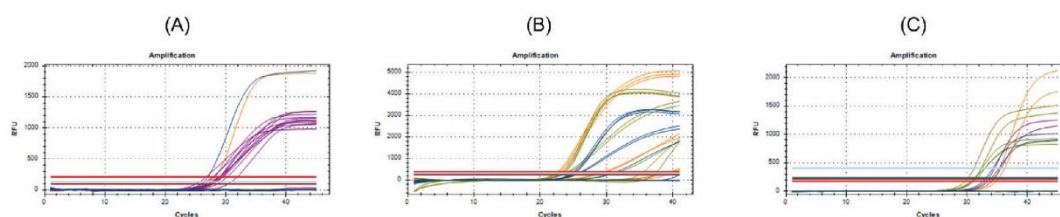
Figure 1. qRT-PCR compatibility testing for COVID-19 Detection. Nasopharyngeal swab and throat swab from patients. After that, 5 μ l of template was sequentially added in the qRT-PCR reagent which was prepared by following the instructions from each manufacture.

DirectAmp™ lysis solution is compatible with multiple fluorophores without a background signal and an interference assay

Colorimetric RT-LAMP Compatibility for COVID-19 Detection

DirectAmp™ lysis solution is compatible with various different swab types including nasopharyngeal (NP) swab, nasal swab and oropharyngeal (throat) swab. The buffer does not interfere the reagent color and performance of RT-LAMP reaction.

Figure 2. Patient sample was collected from (B) nasopharyngeal swab, (C) nasal swab and (D) throat swab was tested with Fastproof™ 30 min-TTR SARS-CoV-2 RT-LAMP Kit (Zenostic, catalog ZE-P14). 5 μ l of template was added in the RT-LAMP reaction. The assay can detect a viral RNA in sample with different sources after incubated at 65°C for 20 mins. Pink color indicated a negative result, while yellow indicated a positive results in which SARS-CoV-2 RNA was detected. (A) positive control assay and (E) negative control assay by using blank DirectAmp™ lysis solution (no template).

Precautions and Disclaimer

- For Research Use Only. Not for use in diagnostic procedures.
- The sample should be kept in range 4°C to -20°C after heating steps to avoid nucleic acid degradation.
- Follow all appropriate biosafety procedures when handling biological samples.

Storage and Stability

DirectAmp™ lysis solution should be stored at 4°C for 3 months

Ordering information

Category	Product	Package volume	cat. No.
Extraction buffer	DirectAmp™ lysis solution	0.3 mL x 100 tubes	ZE-P15

Contact Us
Zenostic Co., Ltd.
Venture Club Bldg. 272 Rama VI Road, ThungPhayathai,
Ratchathewi, Bangkok 10400, Thailand
Tel: +66-2-201-5855 / Fax: +66-2-201-5842
E-mail : info@zenosticbio.com

©2022 Zenostic Co., Ltd. All Rights Reserved.
DirectAmp is a trademark of Zenostic company limited
Product maybe covered by pending or issued patents or
may have certain limitations. Please visit our
Web site for more information

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